KIT VB RAL 10**BLEU DE METHYLENE PUR > 98,5%**

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Principle:

The purpose of the stain test is to assess the clay content of fines by calculating the methylene blue value.

This kit consists of ready-to-use doses for preparing methylene blue solutions at the 10g /L concentration required by European Standard EN 933-9. Each dose has a net weight equivalent to 10 g of dry powder, i.e a weight of $[(100 + W)/10] \pm 0.01$ g where W represents the water content.

Kit description:

- 6 doses equivalent to 10 g of dry powder for preparing 6 solutions of 1 l each at 10 g/liter
- 6 spatulas
- 100 filters with a diameter of 110 mm meeting the standards in force

Specific material required but not provided:

one liter beaker; mechanical stirring device with heating (one liter capacity); one liter volumetric flask

Use:

For professional use only.

Storage temperature: 15-25°C away from light.

Shelf life opened bottle: the solution should not be used more than 28 days after prepared.

Recommendations and/or notes of use:

Follow the method of Preparation of 10 g/liter Methylene Blue Solution in steps C.1.4 to C.2 of annex C of European Standard EN 933-9, i.e.

Bibliography:

British Standard EN 933-9, Assessment of fines. Methylene Blue test.

Standard NF P-94-08, Measuring of the methylene blue adsorption capacity of a rocky soil.

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VB RAL 10 KIT**METHYLENE BLUE MINIMUM PURITY 98.5%**

FOR ASSESSING THE METHYLENE BLUE VALUE BY MEANS OF THE STAIN TEST

Staining Procedure:

Please read all this information carefully before using this device.

STEP 1: Warm 500 ml to 700 ml of distilled or demineralised water in a beaker to a temperature not exceeding 40°C.

STEP 2: Agitate the content of the beaker whilst slowly pouring the methylene blue powder into the warm water. Be sure to have taken all the powder. It is advised to use the pastula enclosed or even rinse the pot with distilled or demineralised water.

STEP 3: Keep agitating for 45 minutes, until complete dissolution of the powder, then allow to cool down to about 20°C.

STEP 4: Pour into a volumetric flask of 1 l capacity, rinse the beaker with distilled or demineralised water to ensure complete transfert of all stain into the flask. Make sure that the flask and the water are at temperature of 20°C and add distilled or demineralised water up to the 1 l graduation mark.

STEP 5: Pour into a tinted glass conservation bottle. The following shall be marked on the conservation bottle:

- 10 g/l methylene blue solution
- Date of preparation
- Limit date of use.

Results:

If observed results vary from expected results please contact RAL DIAGNOSTICS technical service for assistance.

Photo non contractuelle

Non contractual photo

Methylene blue, pure, > 98,5% - 363135-0010

FR: ATTENTION.
EN: WARNING.
DE: ACHTUNG.
ES: ATENCION.
IT: ATENZIONE.
PT: ATENCAO.
AR: امانت.
BG: ВНИМАНИЕ.
CS: POZOR.
DA: OPMÆRKSOMHED.
EL: ΠΡΟΣΟΧΗ.
ET: TÄHELEPANU.
FI: HUOMIO.
HR: PAŽNJA.
HU: FIGYELEM.
JA: 注目.
KO: 주의.
LT: DĖMESIO.
LV: UZMANĪBU.
NL: ATTENTIE.
NO: Attention.
PL: UWAGA.
RO: ATENȚIE.
RU: ВНИМАНИЕ.
SK: POZORNOST'.
SL: POZOR.
SR: ПАЖИЊА.
SV: VAR UPPMÄRKSAM.
TR: DİKKAT.
UK: УВАГА.
ZH: 注意.